

# **GW6010 - GW7010 - GW7015**

**GW6010**: glassware washers with forced-air drying system. hinged door. **GW7010**: Laboratory glassware washer, with forced-air drying system, sliding doors. **GW7015**: Laboratory glassware washer, with forced-air drying system, sliding doors. [ 90 cm wide ]

## - PROGRAM TABLE - (GW)

#### **WARNING**

This document is an appendix to the appliance's user manual and must be consulted together with the manual for proper, safe use of the device.

It contains up-to-date information concerning the characteristics of the washing programs installed in the appliance.

Programs may vary depending on the firmware version installed in the appliance. If in doubt, have authorised staff check on the appliance that the installed version is the same as the version stated in the table at the back of this document.

#### ABBREVIATIONS USED IN THE TABLE

**Demi, DW**: demineralised water, the type of water used in a specific phase of the program.

**Cold, CW**: cold mains water, the type of water used in a specific phase of the program.

**Hot, HW**: hot mains water, the type of water used in a specific phase of the program.

**K\_TEST:** indicates the predisposition in the phase of the "conductivity check" (active only if present and activated the option for the control of the water conductivity).

**ID** prog: identity number of the washing program; it is not necessarily bonded to the program name. (E.g. the program "custom 1" has ID prog = 21).

**Int.:** "Intensive" – use this program to process dirty elements.

N.C.: "NOT CONTROLLED" - refers to the relatively short heating phases during which the heating elements are "on" but no target temperature value is set.

P1, ... P5: the code refers to the activation of the relative peristaltic pump, with dispensing of the liquid detergent associated to it.

Each pump has a default dispensing rate set in the factory.

P1b, P2a: the code refers to dispensing by pump P1, P2 at a rate different from the default rate, used in special cases, as specified in the program table.

PUMP DOSAGE ID.	DEFAULT DOSAGE LABORATORY PROGRAMS GLASSWARE WASHER GW [ml/litro]
P1	5
P1b	3
P2	3
P2a	4
Р3	10
P4	0.4
P1(20), P2(20), P3(20), P4(20)	20 (dosage for service purpose only – peristaltic pump service program)

**TD**: thermal disinfection - phase during which the target temperature is kept above a target value for a specified time. The efficacy of the thermal disinfection process is indicated by means of parameter **A0**. In preset programs, the target value for a thermal disinfection process is always at least 80°C.

**Temp. T** [°C]: Target temperature of the current phase. The temperature inside the tank is kept constant around the specified value for the time shown in the table. The phase during which the temperature is held at the target value is called the **EXTENSION** phase. When no target temperature is indicated or the indication is ("0"), the heating elements are not activated (e.g. during a cold prewash or for cold rinses).





### **VARIOUS NOTES**

**WATER FILLING**: at each step corresponds a load of water. The quantity are specified in the table.

GW6010, GW7010	GW7015
23 liters	32 liters

**PHASE NUMBERING:** the display of the machine, during the progression of the cycle, provides information on the various stages in progress. The first phase assumes the identity of "Phase 1", the second phase "Phase 2", etc.

**CYCLE TIME**: the stated time (on the display) for the complete cycle is completely indicative and refers to standard test conditions: intake water pressure 3bar, cold water temperature 25°C, electrical connection 3/N/PE 400V. Pressures and temperatures lower impact on cycle duration.

### PARAMETER A<sub>0</sub>

Parameter  $A_0$  is used by product standard EN ISO 15883-1 to assign a numerical value to the thermal disinfection process. It is calculated mathematically using the formula:

$$A_0 = \sum 10^{\left[ \left( T - 80 \right) / z \right]} \times \Delta t$$

 $z = 10^{\circ}C - t = duration of time interval considered in seconds - T = load temperature in °C.$ 

When calculating the parameter, only the time intervals during which the temperature is above 65°C are considered.

For our thermal disinfection programs, the calculation is simplified by only including the "extension" phase, when the temperature is kept constant with reference to the target value set.

**DRYING:** Drying stage at the end of the cycle are present only on the models equipped with a drying system.

**CHAMBER RECOVERY DRAIN**: Only if installed the optional "WD-VDS6010": according to the default factory programs settings, the first two phases of the cycle are marked as "recovery drain" (potentially contaminated).

### **FACTORY AND CUSTOM PROGRAMS**

Programs from n.1 to 20 (ID\_prog from 1 to 20) are defined as Factory non-editable programs.

Custom programs: programs from n. 21 to 40 (ID\_prog from 21 to 40) are predisposed to be personalized, with a maximum of n.10 washing phases.

The change can be made through the touch-screen interface of the device, using the updated firmware.

If parameters are modified from the default settings, a record should be kept of the various phases and the new program settings.

**ATTENTION:** The editing of a washing program requires specific knowledge in relation to the load-treatment process, and to the device parameters. For this reason, the function is protected by password.



Always proceed according to the rules in force in the place of installation: **a custom program**, used to process the load, **must be validated according to the standards and regulations applicable**.

Consult the authorized technical service for clarification. Warning: when creating a custom program, do not include more than two thermal disinfection phases at 93°C for 10min.

### **SERVICE PROGRAMS**

Programs 200, 201, 202, 203 are service programs, not to be used for the processing of the load.

19 592 0349 05 GW6010 – GW7015 PROGRAM TABLE 2





# **LABORATORY – PROGRAM TABLE**

		Prewashin	g / Washing / washing	NAOH		NAOH washin o disinfection	ng /	Neutraliza	ition		Mains	rinsing		Demi rin	sing		Rins. Demi ho	ot	Drying	g phase			
ID_prog	Laboratory programs	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C Extension (min)	Temperature °C	Extension (min)	A0	Cycle time std 6010- 7010 [min]	Cycle time std 7015 [min]
1	1 Prewash GW	Cold	0	5																	-	10	12
2	2 Plastic	Cold	0	3	Cold	75	3	Hot P2	NC	3	Hot	NC	1	Demi	NC	1	Demi [K_TEST]	75 1	80	30	-	90	102
3	3 Glassware - quick				Hot P1	75	3	Hot P2	NC	2				Demi	NC	1	Demi [K_TEST]	75 1	80	30	-	70	78
4	4 Glassware - med.				Hot P1	75	5	Hot P2	NC	2				Demi	NC	1	Demi [K_TEST]	75 1	80	30	-	73	81
5	5 Glassware - std.				Hot P1	80	5	Hot P2	NC	2	Hot	NC	1	Demi	NC	1	Demi [K_TEST]	75 1	90	30	-	80	90
6	6 Glassware - int.	Hot	NC	3	Hot P1	90	3	Hot P2ª	NC	2	Hot	NC	1	Demi	NC	2	Demi [K_TEST]	75 1	100	30	1800	88	100
7	7 Glassware (blood)	Cold P1		3	Cold P1	93	3	Hot P2	NC	3	Hot	NC	1	Demi	NC	2	Demi [K_TEST]	75 1	100	30	3600	96	108
8	8 Glassware AGAR	Hot P1	NC	3	Hot P1	93	3	Hot P2	NC	3	Hot	NC	1	Demi	NC	2	Demi [K_TEST]	75 1	100	30	3600	88	100
9	9 AGAR Intensive	Hot P1 P3	80 P4	3	Hot P1 P3	93 P4	5	Hot P2	NC	2	Hot	NC	1	Demi	NC	2	Demi [K_TEST]	75 1	100	30	6000	96	108
10	10 Thermal disin. GW 93°C	11 13	1 1 1		Hot P1	93	3	Hot P2	NC	2				Demi	NC	2	Demi [K_TEST]	75 1	100	30	3600	80	88
11	11 Thermal disin. GW 93°C				Hot P1	93	10	Hot P2	NC	2				Demi	NC	2	Demi [K_TEST]	75 1	100	30	12000	84	92
12	12 Thermal disin. GW int. 93°C 5'	Cold P1	NC	2	Cold	93	5	Hot P2	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	6000	97	109
13	13 Thermal disin. GW int. 93°C 10'	Cold P1	NC	2	Cold	93	10	Hot P2	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	12000	101	113
14	14 Stained Oil	Hot	NC	2	Hot P1 P3	93 P4	1	Hot P2	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	1200	83	95
15	15 Stained oil - int.	Hot P1 <sup>b</sup>	NC	2	Hot P1 P3	93 P4	5	Hot P2	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	6000	89	101
16	16 Oil grease	Hot	50	2	Hot	93	1	Hot P2	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	1200	94	106
17	17 Mineral Oil	Hot	75	1	P1 P3	P4 93	10	Hot	NC	2	Hot	NC	2	Demi	NC	1	Demi [K_TEST]	75 1	100	30	12000	107	119
		P1 P3	P4		P1 P3	P4		P2															





		Prev		g / Wash washing	ing / NA	ОН			IAOH wa disinfect			Neutr	alization		Mains ri	nsing		Demi rin	sing		Rins. Hot d	lemi		Dryir	ng phase			
ID_prog	Laboratory programs		H2O load		Temp °C	Extension (min)		H2O load		Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C Extension (min)	-	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	Temperature °C	Extension (min)	Α0	Cycle time std 6010-7010 [min]	Cycle time std 7015 [min]
18	18 Petrol							Hot		93 5	5	Hot	NC	2	Hot	NC 2		Demi	NC	2	Demi [K_TEST]	75	1	100	30	6000	82	92
							P1	P3	P4			P2																
19	19 Diesel		Hot		NC	3		Hot		93 1	<u> </u>	Hot	NC	2	Hot	NC 2	$\perp$	Demi	NC	2	Demi [K_TEST]	75	1	100	30	1200	95	107
13	15 Diesei	P1	Р3	P4			P1	Р3	P4			P2														1200		107
20	20 Hairman   Datual		Hot		93	10		Hot		93 1	0	Hot	NC	2	Hot	NC 2		Demi	NC	1	Demi [K_TEST]	75	1	100	30	12000	422	125
20	20 Universal Petrol	P1	Р3	P4			P1	Р3	P4			P2ª														12000	123	135

# **CUSTOM PROGRAMS**

On laboratory glassware washers, all the 20 pre-set custom programs have the same structure (as shown in the table below).


		Prewashing , NAOH w			Washing / NA di	AOH wa: isinfecti		Therm	10	Neutral	lization		Ma	ains rinsin	g	Demi r	insing		Rins. Ho	t demi		Drying	phase	
ID_prog	Custom: Program name	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	Temperature °C	Extension (min)	H2O load	Temp °C	Ao
21	Custom 1	DW (32)	-	1																		-	-	-
22	Custom 2	DW (32)	-	1																		-	-	-
23	Custom 3	DW (32)	-	1																		-	-	-
		DW (32)	-	1																		-	-	-
40	Custom 20	DW (32)	-	1																		-	-	-




### **SERVICE PROGRAMS**

		Prewashing NAOH v			Washing / NAOH wa	ashing /	Therm	o disinfe	ection	Neutral	zation		М	ains rinsing		Demi	rinsing		Rins. Hot	demi		Drying	phase			
ID_prog	Service Programs: Program name	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	H2O load	Temp °C	Extension (min)	Temperature °C	Extension (min)	H2O load	Temp °C	Ao	Cycle time std 6010-7010 [min]	Cycle time std 7015 [min]
200	Empty boiler																					-	-	-	-	-
201	Drying																					100	30	-	30	30
202	Service 1 (peristaltic pump test)				CW P1(20)+P3(20)	-	0			CW P2(20)	-	0				CW P4(20)	-	0	DW	-	1	-	-	-	15	15
203	Probes calibration (TL, TCL)	DW	93	10																		90	10	-	31	33

### Recommended use of the service programs:

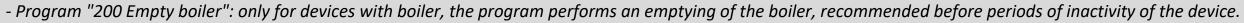

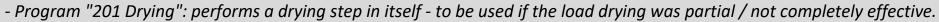



- Program "202 Service 1" It performs a loading of all installed peristaltic pumps can be used to fill the circuit of detergents in case of first installation or to check the proper functioning of the peristaltic (detergents) pumps.
- Program "203 Probes calibration" to verify the good calibration of temperature probes in chamber.

### Please note:

These programs can be disabled by the Superuser to facilitate the choice of the basic operator for the everyday programs.





### **CUSTOM PROGRAM RECORD TEMPLATE**

		Phase 1			Phase 2			Phase 3			Phase 4			Phase 5			Phase 6			Phase 7			Phase 8			Phase 9		l	hase 10	)	Dry	ing	
Name	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Water Filling	T target [°C]	Holding Time [min]	Drying Temperature [°C]	Holding Time [min]	A0 value



Doc.	rev.	Firmware Main	Date	Note
	01	Custom program set cod. 16 813 0000 00	22 / 11 / 2016	2. Custom programs;  3. Service Programs
		Program set cod. 16 812 0013 <b>03</b>		Table structured in three separate sets:  1. Factory (not editable) programs;
19 592 0349 – EN	03	-	20 / 12 / 2016	GW7010.
	04	Program set cod. 16 812 0013 <b>04</b>	18 / 01 / 2017	Ins. Cycle time.
	03	GW7015: Program set cod. 16812002500 Custom program set cod. 16813000400	12/12/2018	GW7015 and conductivity check.
	05	GW6010, GW7010: Program set cod. 16812001305 Custom program set cod. 16813000002	12/12/2019	CW701E and conductivity chack

# Smeg S.p.A. - Instruments Division

Via Leonardo da Vinci, 4 – 42016 Guastalla (RE) Italy Tel. +39 0522 8211 – Fax +39 0522 821 592

E-Mail: <u>instruments@smeg.it</u> - <u>service.instruments@smeg.it</u> <u>www.smeg-instruments.com</u>